


TVFC 400 DOWNLOADABLE DATA LOGGERS PREDOMINANTLY USED BY PROVIDERS

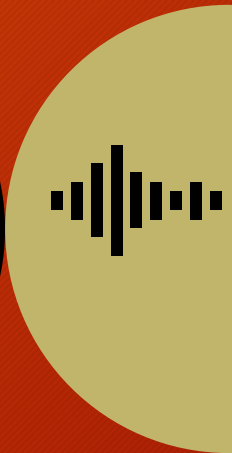

<https://www.vfcdataloggers.com/content/Part%204%20VFC400%20Operation.pdf>

LOGGER COMES WITH GLYCOL BOTTLE

DOWNLOADING CRADLE

ALL DATA LOGGERS THAT MONITOR TVFC/ASN VACCINES MUST HAVE DOWNLOADING CAPABILITIES



CONTROL SOLUTION VFC400 DATA LOGGER KITS

KITS ARE SUPPLIED WITH DATA LOGGER & CRADLE AND CALIBRATED FOR EITHER FREEZER OR REFRIGERATOR - SPECIFIED AT PURCHASE.



DATA LOGGERS WILL HAVE AN EQUIPMENT ID NUMBER LOCATED ON THE BACK, ALWAYS VERIFY EQUIPMENT ID TO THE NUMBER ON THE PROBE'S CORD TO ENSURE THEY READ THE SAME BEFORE CONNECTING THE CORRECT CORD BEFORE POWERING ON.



<https://www.vfcdataloggers.com/vfc400-video-help-guide>

DATALOGGER CERTIFICATE OF CALIBRATION

Certificate of Calibration

Certificate: CSI-175300

Customer Information

Customer: Southwest Pediatric Associates
Address: 7900 FM 1826, Suite 240
City, State Zip: Austin TX 78737

Equipment Information

Equipment ID: V22-13632
Model No: VFC400
Manufacturer: LOG TAG
Serial No: 7862993632
Location: -

Calibration Summary

Calibrated: 11/11/2022
Cal Due Date: 11/11/2024

Temp: 70F +/- 10
Humidity: 50% +/-20
Technician: Lionila De La Cerda Bldg 2

As Found: In Tolerance
Result: Pass

Remarks:



DATA LOGGERS THAT MONITOR TVFC/ASN VACCINES MUST HAVE A CALIBRATION CERTIFICATE - CALIBRATION ENSURES THAT LOGGERS ARE FULLY TESTED AND WORK OPTIMALLY TO CAPTURE ACCURATE TEMPERATURE READINGS.



FOR A CALIBRATION CERTIFICATE TO BE VALID IT MUST HAVE THE FOLLOWING: CUSTOMER INFORMATION, EQUIPMENT INFORMATION, CALIBRATED AND NEXT CALIBRATION DATES.

REMINDER TVFC/ASN ENROLLED CLINICS ARE REQUIRED TO HAVE DATA LOGGERS WITH DOWNLOADING CAPABILITIES FOR ALL UNITS THAT HOLD ANY TVFC/ASN VACCINES.

BACK UP LOGGER REQUIREMENTS

▶▶▶ CLINICS ARE REQUIRED TO HAVE AT LEAST ONE BACK UP DATA LOGGER. IF YOU HAVE MORE THAN ONE REFRIGERATOR AND FREEZER YOU ARE REQUIRED TO HAVE ENOUGH BACKUPS TO REPLACE LOGGERS THAT ARE EITHER UP FOR RECALIBRATION OR ENOUGH LOGGERS TO MONITOR VACCINES DURING EMERGENCY TRANSFERS.

▶▶▶ IT IS BEST TO SEND OUT SOON TO EXPIRE LOGGERS AT LEAST TWO MONTHS BEFORE EXPIRATION. REPLACED DATA LOGGER CERTIFICATES MUST BE SENT TO YOUR R.E. - PLEASE INDICATE ON CERTIFICATE WHAT REFRIGERATOR OR FREEZER AND LOCATION IT IS REPLACING.

▶▶▶ PLEASE ENSURE THAT LOGGER EXPIRATION DATES ARE STAGGERED - MEANING DIFFERENT DATES OF EXPIRY.

NEW OR RE-CALIBRATED DATA LOGGERS

**SUBMIT COPY OF THE CALIBRATION CERTIFICATE TO
YOUR R.E.**

**NOTATE ON CERTIFICATE TYPE OF UNIT
REFRIGERATOR OR FREEZER AND THE LOCATION OF THE
UNIT IN YOUR CLINIC.**

**THE CERTIFICATE WILL BE EMAILED BY R.E. TO DSHS TO
UPDATE STORAGE INFORMATION IN SYNTROPI.**

**VFC DATALOGGER PURCHASING & RE-CALIBRATION
CONTROL SOLUTIONS - <https://.vfcdataloggers.com/>**